

# Local Run Manager v3

## Software Guide

Local Run Manager Overview	3
Off-Instrument Installation	4
User Passwords	7
Dashboard Overview	9
Administrative Settings and Tasks	13
Workflow Overview	24
Analysis Module Updates	29
Revision History	30
Technical Assistance	31



This document and its contents are proprietary to Illumina, Inc. and its affiliates ("Illumina"), and are intended solely for the contractual use of its customer in connection with the use of the product(s) described herein and for no other purpose. This document and its contents shall not be used or distributed for any other purpose and/or otherwise communicated, disclosed, or reproduced in any way whatsoever without the prior written consent of Illumina. Illumina does not convey any license under its patent, trademark, copyright, or common-law rights nor similar rights of any third parties by this document.

The instructions in this document must be strictly and explicitly followed by qualified and properly trained personnel in order to ensure the proper and safe use of the product(s) described herein. All of the contents of this document must be fully read and understood prior to using such product(s).

FAILURE TO COMPLETELY READ AND EXPLICITLY FOLLOW ALL OF THE INSTRUCTIONS CONTAINED HEREIN MAY RESULT IN DAMAGE TO THE PRODUCT(S), INJURY TO PERSONS, INCLUDING TO USERS OR OTHERS, AND DAMAGE TO OTHER PROPERTY, AND WILL VOID ANY WARRANTY APPLICABLE TO THE PRODUCT(S).

ILLUMINA DOES NOT ASSUME ANY LIABILITY ARISING OUT OF THE IMPROPER USE OF THE PRODUCT(S) DESCRIBED HEREIN (INCLUDING PARTS THEREOF OR SOFTWARE).

© 2022 Illumina, Inc. All rights reserved.

All trademarks are the property of Illumina, Inc. or their respective owners. For specific trademark information, see [www.illumina.com/company/legal.html](http://www.illumina.com/company/legal.html).

## Local Run Manager Overview

The Local Run Manager software is an integrated solution for recording samples for a run, specifying run parameters, monitoring status, analyzing sequencing data, and viewing results.

Local Run Manager has the following features:

- ▶ Integrates with the instrument control software when run on the instrument computer.
- ▶ Records samples and input files, such as manifests and reference genomes.
- ▶ Provides a run setup interface specific to the selected analysis module.
- ▶ Performs data analysis on completed sequencing runs.
- ▶ Performs a series of analysis steps specific to the selected analysis module.
- ▶ Shows analysis metrics upon completion of analysis.

Only a single instance of Local Run Manager can be installed on the computer.

## Analysis Modules

The Local Run Manager software uses different analysis modules to perform data analysis. Each module has a specific set of procedures appropriate for an application and library type. For a description of each analysis module, see the [Local Run Manager support page](#) on the Illumina® website.

The Local Run Manager software does not include any preinstalled analysis modules. Each analysis module, and the off-instrument version of Local Run Manager, must be downloaded and installed on your system separate from Local Run Manager. Not all modules can be used with the off-instrument version of Local Run Manager. For more information on installing an analysis module, see the Local Run Manager Workflow Guide specific to that module.

## Viewing Local Run Manager

The Local Run Manager interface is viewed through a web browser. Open a web browser on a computer with access to the same network as the instrument and connect using either of the following methods:

- ▶ On a local network, connect using the instrument name. For example, **https://myinstrument**.
- ▶ To view the Local Run Manager interface on the instrument monitor, open a web browser and enter **https://localhost** in the address bar.
- ▶ To best view the Local Run Manager interface on a computer separate from the instrument, make sure that the zoom setting on your browser is set to 100%.
- ▶ To access the instrument remotely, port 80 must be **open**.

## Sequencing During Analysis

The instrument computing resources are dedicated to either sequencing or analysis.

If a new sequencing run is started before secondary analysis of an earlier run is complete, a confirmation dialog box appears. After confirming you want the new sequencing run to start, Local Run Manager stops secondary analysis of the earlier run until the new run completes sequencing.

After the new run completes sequencing, secondary analysis of the earlier run automatically starts again from the beginning.

## Off-Instrument Installation

The instructions included in this section only apply to the off-instrument installation of Local Run Manager. For software issues on your Illumina instrument, refer to the system guide for your instrument.

### Requirements

To use the *off-instrument* implementation of Local Run Manager, your computer must meet the following requirements.

#### Computing Requirements

- ▶ Minimum 16 GB RAM
- ▶ Minimum 1 TB disk space
- ▶ 64-bit quad core processor (2.8 GHz or higher) with the AVX instruction set (recommended) or Intel Core i7-210QE 2.10 GHz (or equivalent) processor with the AVX instruction set (minimum)

#### Software Requirements

- ▶ Windows 10 (Personal Edition version is not supported)—For more information, see [Change System Language Settings](#) on page 6
- ▶ Microsoft .NET 4.5.2 or later
- ▶ Chrome 76.0.3809.132-x64 or later
- ▶ Internet Information Services

### Dependencies Installed by Local Run Manager

Uninstall all components from the previous version of Local Run Manager before proceeding to install version 3 of Local Run Manager. Failure to do so will cause a delay and a complaint will need to be initiated.

If any of the following components are already present on your system, installation may fail.

- ▶ PostgreSQL 9.3
- ▶ Rabbitmq 3.5.1
- ▶ Erlang/OTP 17
- ▶ C++ 2012 64-bit Redistributable
- ▶ C++ 2013 64-bit Redistributable
- ▶ C++ 2015 64-bit Redistributable
- ▶ Illumina Run Copy Service (applicable to off-instrument installations)
- ▶ Internet Information Services
- ▶ Local Run Manager

### Install Off-Instrument Local Run Manager

- 1 Uninstall the Local Run Manager v2.
- 2 Navigate to Programs & Features.
  - ▶ Ensure the Erlang, Postgres, RabbitMQ, and Local Run Manager are uninstalled.

- 3 Navigate to %AppData% in Windows Explorer.
  - ▶ Delete the RabbitMQ folder under Roaming folder. If you are unable to delete it, open Task Manager, end the task for epdm.exe, and restart the PC.
- 4 Navigate to C:\Program Files.
  - ▶ Delete the PostgreSQL folder.
  - ▶ Delete erl6.4 folder. If not able to delete it, open Task Manager and end the task for erl.exe and ertsrv.exe, and restart the PC if necessary.
- 5 Navigate to C:\Program Files (x86)
  - ▶ Delete the RabbitMQ Server folder.
- 6 Launch Services.
  - ▶ Stop Illumina Local Run Manager Analysis Service.
  - ▶ Stop Illumina Local Run Manager Job Service.
- 7 Launch the Registry Editor.
  - ▶ Navigate to HKEY\_LOCAL\_MACHINE\SYSTEM\CurrentControlSet\services.
    - ▶ Delete RabbitMQ.
    - ▶ Delete LocalRunManagerAnalysisService.
    - ▶ Delete LocalRunManagerJobService.
    - ▶ Delete LocalRunManagerImpersonationService.
  - ▶ Navigate to HKEY\_LOCAL\_MACHINE\SOFTWARE\Ericsson\
    - ▶ Delete Erlang.
  - ▶ Navigate to HKEY\_LOCAL\_MACHINE\SOFTWARE\WOW6432Node\Ericsson
    - ▶ Delete Erlang.
- 8 Navigate to C:\Users\<Windows user account that installed RabbitMQ>\
  - ▶ Delete .erlang.cookie.
- 9 Navigate to C:\Windows\
  - ▶ Delete .erlang.cookie.
- 10 Launch Environment Variables. Under System variables,
  - ▶ Delete ERLANG\_HOME
  - ▶ Delete RABBITMQ\_NODENAME
- 11 Remove or rename the following folders:
  - ▶ C:\Illumina\Local Run Manager
  - ▶ C:\Illumina\Local Run Manager Analysis Job Service
  - ▶ C:\Illumina\Local Run Manager Analysis Service
  - ▶ C:\Program Files\Illumina\Local Run Manager
  - ▶ C:\ProgramData\Illumina\Local Run Manager Analysis Job Service
  - ▶ C:\Program Files\Illumina\Local Run Manager Analysis Service
  - ▶ C:\ProgramData\Illumina
  - ▶ C:\ProgramData\Illumina\Local Run Manager Analysis Service
  - ▶ C:\ProgramData\Package Cache
- 12 Delete localrunmanager registries in Registry Editor.
- 13 Right-click **OffInstrument\_Research\_(version).exe** **Run as administrator** and select **Run as administrator**.
- 14 Select **Install LRM v3**.

15 When installation is complete, select **Close**.

## **Change System Language Settings**

To use Local Run Manager, the system language must be set to English (United States). To change the system language settings for your system, see the reference guide for your operating system.

## User Passwords

The user management option, which requires password verification to use Local Run Manager, is disabled by default. But if your system is secured with this option enabled, you need a valid user name and password to log in to the system. Only an admin user can assign user credentials.



### NOTE

User accounts are specific to an instrument. User passwords are not universal across multiple instruments.

A message appears at the top of the interface to remind you to reset your password when it has expired.

## My Account

You can reset your existing password at any time using the My Account window. From My Account, you can view your user name, assigned role, and permissions. You can also configure your security questions to regain access to your account in case you get locked out of Local Run Manager.

To change your password, you must know your current password. If you forget your password, you can answer your account's security questions. If you cannot answer the security questions correctly, an Admin can assign a temporary password so you can update your password.

## Password States

The following are possible password states:

- ▶ **Temporary password**—When an Admin creates a user account, the Admin assigns a temporary password for the new user.
- ▶ **User password**—Upon first access, a new user is prompted to change the temporary password from the login screen to a user-assigned password. The user must also set up security questions to help verify that the user is the person requesting access to the account.
- ▶ **Forgotten password**—If a user forgets a password, an Admin can assign a temporary password that can be changed upon next access.
- ▶ **Used password**—Users cannot reuse a password for up to five password cycles.
- ▶ **User lockout**—An Admin can change the limit for number of attempts to log in using an invalid password. The default is 5. If a user exceeds the number of allowed attempts, the user account is locked. Only an Admin can unlock the account, or assign a temporary password.

## Reset Your Password

Follow the steps below to reset a password.

- 1 From the navigation bar, select the **User** icon next to your name.
- 2 From the drop-down list, select **My Account**.
- 3 In the Password section, select **Edit**.
- 4 In the Old Password field, enter your old password.
- 5 In the New Password field, enter a new password.
- 6 In the Confirm New Password field, reenter your new password.
- 7 Select **Save**.

## Disable Password Expiration

Disabling password expiration is only available to admin users. Follow the steps below to disable expiration dates for passwords.

- 1 From the navigation bar, select the Tools drop-down list and then select **System Settings**.
- 2 Select the Security tab.
- 3 Under Password Expiration, clear the **Disable password expiration** checkbox.

## Set Up Security Questions

When logging into Local Run Manager for the first time, users are asked to set up security questions. If a user forgets their password, correctly answering their account's security questions will unlock the account. Follow the steps below to set up security questions for the account.

- 1 From the first Question drop-down menu, select a security question. You can also make a custom question by selecting **Create my own question** and entering a question of your own.
- 2 Enter the answer in the Answer field.  
Answers are not case-sensitive.
- 3 Repeat steps 1 and 2 for the remaining question and answer fields.
- 4 Select **Update Security Questions** to save the security information to your account.

Security question information can be changed later by going to the My Account menu and selecting **Reset** in the Security Questions section.



## Dashboard Overview

When you log in to the Local Run Manager software, the dashboard opens. From the dashboard, you can perform the following tasks:

- ▶ Track sequencing and analysis status
- ▶ Sort and filter runs
- ▶ Create and edit runs
- ▶ Hide runs
- ▶ Import runs
- ▶ Stop or requeue analysis
- ▶ View analysis status
- ▶ Manage run folder

The dashboard lists all runs performed on the instrument, 10 runs per page. Use the navigation arrows at the bottom of the list to scroll through pages.

Each run listed includes the run name, analysis module, run status, and the date that the run was last modified.

- ▶ **Run Name/ID**—Lists the assigned run name and links to the run results pages. The Run ID is added to the run name column at the start of sequencing or analysis.
- ▶ **Module**—Lists the analysis module assigned to the run.
- ▶ **Status**—Lists the status of the run and includes a progress status bar. For more information, see *Possible Run States on page 28*.
- ▶ **Last Modified**—Lists the date and time that a run or analysis session was last modified. By default, the Active Runs page is sorted by this column.

To sort the run list, select any column heading to sort in ascending or descending order.

## Active Runs Page

The dashboard opens on the Active Runs page by default. At the top of the Active Runs page, a summary of active runs provides a count of runs in each of the following run status categories.

- ▶ **Ready**—Indicates that the run is ready for sequencing or importing.
- ▶ **In Progress**—Indicates that the run is in the process of sequencing or data analysis, or in the process of stopping after a manual stop command.
- ▶ **Stopped or Unsuccessful**—Indicates that analysis was stopped manually or that analysis was unsuccessful.
- ▶ **Complete**—Indicates that analysis was successfully completed.
- ▶ **Total**—The total number of runs on the Active Runs page.

## Filter Runs

- 1 Select a run status category in the summary of active runs to filter the Active Runs page as follows.
  - ▶ **Ready**—Filters the list to show only runs that are ready for sequencing.
  - ▶ **In Progress**—Filters the list to show only runs that are in progress.
  - ▶ **Stopped or Unsuccessful**—Filters the list to show only runs that were stopped or unsuccessful.
  - ▶ **Complete**—Filters the list to show only runs that are complete.

- ▶ **Total**—Removes filters and shows all active runs.

## Hide and Restore Runs

- 1 From the Active Runs page, select **Actions** next to the run name that you want to hide.
- 2 Select **Hide**.
- 3 When prompted to confirm the move, select **Hide**.
- 4 To show the Hidden Runs page, select the Active Runs drop-down arrow and select **Hidden Runs**.
- 5 To restore a run to the Active Runs page, select **Actions** next to the run name that you want to restore.
- 6 Select **Restore**.
- 7 When prompted to confirm the move, select **Restore**.
- 8 To show the Active Runs page, select the Hidden Runs drop-down arrow and select **Active Runs**.

## Delete a Run Folder

You can manually delete run folders to manage storage space. To delete a run folder, the run must be in one of the following states:

- ▶ Sequencing Complete
- ▶ Primary Analysis Complete
- ▶ Primary Analysis Unsuccessful
- ▶ Sequencing Errored
- ▶ Analysis Errored
- ▶ Analysis Complete
- ▶ Sequencing Stopped
- ▶ Analysis Stopped

- 1 From the Active Runs page, select **Actions** next to the run name that you want to delete.
- 2 Select **Delete Run Folder**.  
The **Delete Run Folder** option is replaced with **Relink**. A trash bin icon appears on the run.

## Relink a Run Folder

Only Administrators can relink deleted run folders.

- 1 Copy the run folder back to its original location.  
If the run is not saved in its original location, the software displays an error message following an attempt to restore.
- 2 Select **Actions** next to the run.
- 3 Select **Relink Run Folder**.
- 4 Select **Confirm** to proceed and relink the run.  
The run is restored to the state it was in before the run folder was deleted.

## Search for Runs or Samples

- 1 From the navigation bar, select the Search icon.  
Searches can only be performed on active runs.
- 2 In the search field, enter a run name or sample ID.

As you type, a list of possible matches appears to assist your search.

- 3 Select a match from the list or press **Enter**.
    - ▶ If you entered a run name in the search field, the Run Overview tab opens.
    - ▶ If you entered a sample ID in the search field, the Samples & Results tab opens.
- For more information, see [View Analysis Results Overview on page 26](#).

## Edit a Run

- 1 From the Active Runs page, select **Actions** next to the run name that you want to edit.
- 2 Select **Edit**.
- 3 When prompted to confirm the action, select **Continue**.
- 4 Edit run parameters as needed.
- 5 When finished, select **Save Run**.  
The Last Modified date for the run is updated on the Active Runs page.

## Import a Run

- 1 From the Active Runs page, select **Create Run**.
- 2 Set the parameters for the run. For more information, see the workflow guide for the analysis module.
- 3 Select **Save Run**.
- 4 From the Active Runs page, select **Actions** next to the newly created run.
- 5 Select **Import**.
- 6 Enter the run folder path of the run that you want to import.  
Runs imported from network locations may take longer to analyze than runs imported from a local drive.
- 7 Select **Import Run**.  
Analysis starts immediately.

## Stop Analysis

- 1 From the Active Runs page, select **Actions** next to the run name that you want to stop.
- 2 Select **Stop Analysis**.  
The Stop Analysis command appears for runs that are either in the state of Analysis Queued or Analysis Running.
- 3 When prompted to confirm, select **Stop Analysis**.

## Requeue Analysis

You might requeue analysis if analysis was stopped, if analysis was unsuccessful, or if you want to reanalyze a run with different settings. You can requeue analysis from the dashboard or from the analysis results screen.

### Requeue Analysis from Active Runs

- 1 From the Active Runs page, select **Actions** next to the run name that you want to requeue.
- 2 Select **Requeue**.

- 3 When prompted, select from the following options:
  - ▶ Select **Edit Setup** to change analysis parameters. From the Requeue Analysis page, change preferred settings, and then select **Requeue Analysis**.
  - ▶ Select **Requeue**. Analysis begins using current analysis parameters.

## Requeue Analysis from Results Page

- 1 Select the run name on the Active Runs page.
- 2 Select **Requeue Analysis**.
- 3 When prompted, select from the following options:
  - ▶ Select **Edit Setup** to change analysis parameters. From the Requeue Analysis page, change preferred settings, and then select **Requeue Analysis**.
  - ▶ Select **Requeue**. Analysis begins using current analysis parameters.

## Change the Output Run Path

You can change the location of the output run folder when a run finishes. This is useful when you want to requeue a run and save the output to a new folder.

- 1 From the Run Dashboard, filter the run list to show the completed runs by selecting **Completed**.
- 2 Select the run that you want to edit.  
Make sure that the selected run is not in progress or the output run folder cannot be changed.
- 3 Select the Run Overview tab.
- 4 From the Output Run Folder section, select **Edit**.
- 5 Enter a new path in the Output Run Folder field and select **Save**.  
You can only edit the path to the output folder, the output folder name remains the same.

## View On-Screen Help

Each page provides help to guide you through the available options on the screen. From the navigation bar, select **Help** for information on the possible choices.

## Administrative Settings and Tasks

The following settings and tasks are available from the Tools drop-down menu:

Tool	Description	Role Permission
Audit Trails	Monitors access frequency and user activity.	Admin
Index & Library Prep Kits	Allows the addition of custom indexes and library prep kits. For more information, see <a href="#">Index &amp; Library Prep Kits on page 21</a> .	User*
Modules & Manifests	Allows upload of manifest files for use with all runs using a specific analysis module.	User*
System Settings	Sets system security parameters and maintenance intervals.	Admin
Account Management	Adds users, assigns permissions, and sets passwords.	Admin

\* Available on a per user basis

## Account Management

All user accounts are listed on the Account Management page, which includes the first name, last name, and user name for each account. This setting is turned off by default.



### NOTE

User accounts are specific to an instrument. User passwords are not universal across multiple instruments.

Within the details of each account, the role and associated permissions are managed. Two roles are possible, either Admin or User.

- ▶ **Admin**—Admin roles have full permissions, by default.
- ▶ **User**—User roles are configurable based on a subset of possible permissions. All users can create runs regardless of permission settings.

Only an Admin can view the Account Management page.

## User Permissions

The following table lists the permissions and roles that are allowed to configure each permission.

Permission	Description	Admin Role	User Role
Manage User Accounts	Create and edit user accounts.	Allowed	Not allowed
Edit Local Run Manager System Settings	Set security and maintenance settings.	Allowed	Not allowed
Upload Site Certificate	Add custom HTTPS certificate for LRM.	Allowed	Not allowed
Relink Run Folder	Restore deleted run data back to a run.	Allowed	Not allowed
View Audit Trails	View, filter, and export audit trails.	Allowed	Not allowed
Requeue Analysis	Reanalyze runs; edit analysis parameters.	Allowed	Optional
Edit Module Settings	Upload manifest files for analysis modules.	Allowed	Optional
Manage Kits	Create a custom library kit definition.	Allowed	Optional

## Create a New User

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **Account Management**.
- 2 From the Account Management page, select **Create User**.
- 3 In the Create New User dialog box, enter the first and last name of the new user.
- 4 In the User Name field, enter a user name.  
User names must be unique and cannot be reused or changed.
- 5 In the Password field, enter a temporary password.  
Temporary passwords are not stored in the password history and can be reused.
- 6 In the Confirm Password field, reenter the temporary password.
- 7 If you are creating this account for someone else, you must set up security questions for the account.  
For more information, see [Set Up Security Questions](#) on page 8.
- 8 To select a role, select **Admin** or **User**.
- 9 Select user permissions based on the specified user role.
- 10 Select **Create User**.

## Reset User Password

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **Account Management**.
- 2 Locate the user name you want to edit, and select **Edit**.
- 3 In the New Password field, enter a temporary password.  
Temporary passwords are not stored in the password history and can be reused.
- 4 In the Confirm Password field, reenter the temporary password.
- 5 Select **Update User**.

## Unlock User

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **Account Management**.
- 2 Locate the user name you want to edit, and select **Edit**.
- 3 From the Edit User dialog box, select **Unlock User**.
- 4 In the New Password field, enter a temporary password.
- 5 In the Confirm Password field, reenter the temporary password.
- 6 Select **Update User**.

## Change User Permissions

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **Account Management**.
- 2 Locate the user name you want to edit, and select **Edit**.

- 3 To change a role, select **Admin** or **User** to toggle between the options.
- 4 Select user permissions based on the specified user role.
- 5 Select **Update User**.

## Disable a User

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **Account Management**.
- 2 Locate the user name you want to edit, and select **Edit**.
- 3 In the Edit User dialog box, select **Disable User**.
- 4 When prompted to disable the user, select **Disable**.

## System Settings

System settings are global parameters for user security, automated data maintenance, service accounts, reference genome repository assignment, and program certificates. The following tasks can be managed in the system settings:

- ▶ System security settings include duration for password expiration, maximum login attempts, and duration for idle timeout. Available when User Management is enabled.
- ▶ System maintenance includes automatic cleanup of inactive run folders, database backup frequency, and a command for immediate database backup.
- ▶ Service account settings include changing user accounts for data analysis, performing backups, and data storage.
- ▶ Assign the reference genome repository location.
- ▶ Manage and download Local Run Manager certificates.
- ▶ Change the network output path.

Most of these settings can only be seen by an Admin but all users are able to download the server certificate to access Local Run Manager with HTTPS.

## Specify System Security Settings

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **System Settings**.
- 2 Select the Security tab.
- 3 From the Password Expiration section, specify the number of days before a password expires and has to be reset or select the **Disable password expiration** checkbox.
- 4 Specify the number of days that a user is reminded to reset a password before it expires.
- 5 From the User Lockout section, specify the maximum number of attempts that the user can enter invalid login credentials before being locked out of the system.
- 6 From the Auto Logout section, specify the number of minutes that the system can remain idle before the user is automatically logged out.
- 7 Select **Save**.

## Specify System Maintenance Settings

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **System Settings**.
- 2 Select the Maintenance tab.
- 3 **[Optional]** Select the **Enable Automatic Deletion** checkbox to enable automatic removal of older analysis files.  
When enabled, the system keeps recent analysis information and deletes all other temp files to improve performance.
- 4 If the Enable Automatic Deletion checkbox is selected, specify the automatic deletion trigger. This parameter varies by instrument and is specified as one of the following:
  - ▶ Folder inactivity (in days)
  - ▶ Folders to keep
- 5 Enter a preferred path to a backup location other than on the local instrument computer, such as a network location or on another networked computer.
- 6 In the Backup Period field, enter the duration in days between each backup.
- 7 To create an immediate backup, select **Backup Now**.
- 8 Select **Save**.

## Specify Service Account Settings

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **System Settings**.
- 2 Select the Service Accounts tab.
- 3 To enable and set the credentials for Analysis Service, Job Service, and Universal Copy Service, select Windows Account.  
The credentials for the account with access to network drive must be saved in Windows Credential Manager



### NOTE

Make sure that the service account you choose has permission to read and write to the output run folder.

- 4 In the User Name field, enter a user name.  
Include .\ at the beginning of the user name (.\username). For domain users, include the domain name, then a backslash in the user name (domain\username).
- 5 In the Password field, enter a password.  
For instruments that run Windows 10, the operating system requires a Windows password change every 180 days. The Windows Account in Local Run Manager must be updated with the same password as the Windows operating system.
- 6 Select **Save**.

## Specify Reference Genome Repository Path

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **System Settings**.
- 2 Select the System Configuration tab.
- 3 In the Repository Path field, enter the path of the reference genome repository.



The repository must be on a local or network folder. Access to online repositories is not supported. You can also select **Set to Default** to reset the path to C:\Illumina\Genomes.

- 4 Select **Save**.

## Set Application Language

Local Run Manager comes installed with the English language pack only. To add other languages, contact Illumina technical support. To set the language for Local Run Manager, perform the following.

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **System Settings**.
- 2 Select the System Configuration tab.
- 3 From the language drop-down list, select the desired language.
- 4 Select **Save**.

## Specify Certificates

Local Run Manager uses HTTPS to encrypt data connections to make sure that run data information is private and secure. Security certificate settings should only be changed in the following situations:

- ▶ Access Local Run Manager from a new remote device
- ▶ Use your own custom security certificate
- ▶ Regain access to Local Run Manager after changing the system name or moving the system to a new domain

## Install a Security Certificate For Remote Access

- 1 On the remote device, login to Local Run Manager.
- 2 From the dashboard navigation bar, select the Tools drop-down list, and then select **System Settings**.
- 3 Select the Certificates tab.
- 4 Select **Download** to download a zip file that contains the Local Run Manager certificate.
- 5 Extract the contents of the zip file.
- 6 Right-click the BAT file and select **Run as administrator**.
- 7 Follow the prompts to finish the installation, and then restart your browser.

## Upload a Custom Security Certificate

To replace the existing Local Run Manager certificate, do the following.

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **System Settings**.
- 2 Select the Certificates tab.
- 3 Select **Browse**, select the new certificate file (\*.pfx), and then select **Open**.
- 4 Enter the password for the certificate file, and then select **Set Certificate**.
- 5 Restart your browser.

## Regenerate a Security Certificate

If there was recent change to the system name or the system was moved to a new domain, you must regenerate the security certificate to regain access to Local Run Manager.

- 1 Visit <https://localhost> in the system browser.
- 2 Enter your user credentials if prompted.
- 3 From the dashboard navigation bar, select the Tools drop-down list, and then select **System Settings**.
- 4 Select the Certificates tab.
- 5 Select **Regenerate the Certificate**.
- 6 Restart your browser.
- 7 If you need to connect to Local Run Manager from another device, download and install the regenerated certificate on the remote device.

## Change the Network Output Path

For some types of instruments, you access the network output path to the output folder through the instrument control software. In other types of instruments, you access the path through Local Run Manager.

In Local Run Manager, navigate to the network output path. Instructions can vary.

- 1 Select **Manage Instrument**.
- 2 Select **System Configuration**.
- 3 Select **Analysis Configuration**.
- 4 Select the **Local Run Manager** tab.  
The default user is admin.
- 5 In the Output folder, enter the full UNC file path to the output folder.  
Make sure to follow UNC path instructions in *Designate the Universal Naming Convention (UNC) Network Output Path* on page 18.
- 6 Select **Save**.

Both the Windows Account in Local Run Manager and the Windows account on the operating system of the instrument require permission to read and write to the output folder. To check permissions, consult your IT administrator. To set up the Windows Account in Local Run Manager, see *Specify Service Account Settings* on page 16.

## Designate the Universal Naming Convention (UNC) Network Output Path

In the output folder, enter the full UNC file path. The UNC path includes two backslashes, the server name and the directory name, but *not* a letter for a mapped network drive.

- ▶ Paths to the output folder that are one level require a trailing backslash.
  - ▶ Example UNC path: \\servername\directory1\
- ▶ Paths to the output folder that are two or more levels do not require a trailing backslash.
  - ▶ Example UNC path: \\servername\directory1\directory2
- ▶ Paths to a mapped network drive cause errors. *Do not use*.
  - ▶ Example of a mapped network drive path: T:\sbsfiles

## Modules & Manifests

The Modules & Manifests page provides a list of installed analysis modules in the left navigation panel. Selecting an analysis module name opens a page that lists the module version and last modified date.

When a new or updated analysis module is available, you can download the module from the Illumina website and install it on the Local Run Manager system. For more information, see [Analysis Module Updates on page 29](#).

For analysis modules that require manifests, use the Add Manifests feature to add manifests to the module. When you add manifests for an analysis module from the Modules & Manifests page, the manifests are available for all runs using that analysis module.

Alternatively, you can import a manifest for a specific run using **Import Manifests** on the Create Run page. When you import a manifest from the Create Run page, the manifest is available for the current run only.

### Add Manifest Files to an Analysis Module

- 1 Make sure that the manifest you want to copy to the module are available in an accessible network location or on a USB drive.
- 2 From the dashboard navigation bar, select the Tools drop-down list, and then select **Modules & Manifests**.
- 3 From the list of analysis modules on the left panel, select the appropriate module name.
- 4 Select **Add Manifests**.
- 5 Navigate to the manifest and select the manifest that you want to add.  
You can add multiple manifests.

### Export a Manifest File

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **Modules & Manifests**.
- 2 From the list of analysis modules on the left panel, select the appropriate module name.
- 3 Select **Export** next to the manifest file that you want to export.

### Delete a Manifest File

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **Modules & Manifests**.
- 2 From the list of analysis modules on the left panel, select the appropriate module name.
- 3 Select **Delete** next to the manifest file that you want to remove.

## Audit Trails

Audit trails records information about specific actions, such as user access, user profile changes, and changes to system, run, or analysis parameters. Each audit trail entry includes the following information:

- ▶ When, represented as date YYYY-MM-DD and time in 24-hour time format.
- ▶ Who, represented by the user name who initiated the action.
- ▶ What, represented by a short predefined description of the action taken.
- ▶ Affected Item, represented by 4 affected item categories of User, Run, Analysis, or System.

To sort the audit trails list, select any column heading to sort in ascending or descending order. You can filter the audit trails list to show specific data and then export the filtered list to a report in PDF file format. Only an admin user can view the Audit Trails page.

## Filter Audit Trails

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **Audit Trails**.
- 2 Select **Filter**.  
You can apply filters to previously filtered results to refine the list even further. To apply a filter to the entire audit trails database, clear previous filters before proceeding.

### Filter by Date

- 1 In the From field, enter a date (in yyyy-mm-dd format) to specify the beginning date for the filtered results. Alternatively, select the **Calendar** icon, and then select a From date.
- 2 In the To field, enter a date (in yyyy-mm-dd format) to specify the end date for the filtered results. Alternatively, select the **Calendar** icon, and then select a To date.
- 3 Select **Filter**.

### Filter by User Name

- 1 In the Who field, enter a user name.  
You can enter any portion of a user name. Asterisks (\*) are not required.
- 2 Select **Filter**.

### Filter by Action

- 1 In the What field, enter an action description.  
You can enter any portion of an action description. Asterisks (\*) are not required.
- 2 Select **Filter**.

### Filter by Affected Item Description

- 1 In the Affected Item field, enter any portion of the affected item description.  
The description can be a run name, user name, analysis module name, or report name, for example. You can enter any portion of a description. Asterisks (\*) are not required.
- 2 Select **Filter**.

### Filter by Affected Item Category

- 1 To filter by an affected item category, select the **Affected Item** icon and select from the following options:
  - ▶ **Analysis**—Filters the list to show changes in analysis parameters or analysis status.
  - ▶ **Run**—Filters the list to show changes to run parameters or run status
  - ▶ **System**—Filters the list to show actions at the system level, such as file uploads, system maintenance, or security settings.
  - ▶ **User**—Filters the list to show user actions and user authentication actions.
- 2 Select **Filter**.

## Export Audit Trails

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **Audit Trails**.
- 2 Select **Filter**.
- 3 Apply preferred filters.
- 4 Select the **Export**.  
The software generates a PDF report that includes the user name, date of export, and filter parameters.

## Index & Library Prep Kits

The Index & Library Prep Kits page provides a list of compatible index and library prep kit definitions installed with the software. A kit description and read type are listed for each kit.

From the Index & Library Prep Kits page, you can upload custom index or library prep kit definition files. Library prep kit definition files list information about the prep kit such as compatible analysis modules, index sequences, and the maximum number cycles or reads in a run. Index kit definition files specify details about the index kit such as index strategy, index sequences, and supported library prep kits. Select **Download Template** from the Library Prep Kit or Index Kit tabs or visit the Local Run Manager support page on the Illumina website for examples of kit definition files.

Guidelines for adding a custom library kit:

- ▶ Adding a custom library prep kit is a feature for advanced users.
- ▶ Always assign a *unique* kit name. Custom library prep kit names cannot be reused on the same instrument.

## Create a Custom Library Prep Kit Definition File

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **Index & Library Prep Kits**.
- 2 From the Library Prep Kit tab, select **Download Template**.  
A library prep kit definition template file, local-run-manager-library-prep-kit-template.tsv, is saved to your computer.
- 3 Open the template file in a text editor and edit the file with your custom parameters.  
The template file has the following section titles, headings, and information. Make sure that you separate the parameters with tabs.
  - ▶ **[Kit]**—In this section, specify a kit name and default run parameters.

Row Heading	Description
Name	Assign a unique kit name for the new kit.
Description	A brief description or long kit name.
IndexStrategy	Options are <b>NoIndex</b> , <b>SingleOnly</b> , <b>DualOnly</b> , <b>SingleAndDual</b> , <b>NoAndSingle</b> , <b>NoAndDual</b> , or <b>All</b> . Enter the index strategy exactly as shown with no spaces.
ReadType	Options are <b>Single</b> , <b>Paired</b> , or <b>All</b> .
DefaultReadLength1	The default number of cycles for Read 1, such as <b>151</b> .
DefaultReadLength2	The default number of cycles for Read 2, such as <b>151</b> .

- ▶ **[Resources]**—In this section, specify custom parameters, such as the maximum number of samples possible.

Column Heading	Description
Name	Enter a name for the parameter. For example, Sample Limit. Spaces are allowed.
Type	Enter a name for the parameter without spaces. For example, SampleLimit.
Format	Enter INT.
Value	Enter a value for the parameter. For example, 384.

- ▶ **[Indices]**—In this section, specify which index sequences are used with the kit.

Column Heading	Description
Name	List the index sequence name, such as <b>A701</b> .
Sequence	List the index sequence.
IndexReadNumber	Specify the Index Read, either <b>1</b> or <b>2</b> .



#### NOTE

During analysis, the iSeq™ 100, MiniSeq™, and NextSeq™ Systems automatically reverse complement the i5 indexes in custom library prep kits. Make sure that the i5 indexes are in the forward orientation.



#### NOTE

Sample sheets produced in Illumina Experience Manager (IEM) are not compatible for use with Local Run Manager. Sample sheets produced in IEM are not compatible for use with Sample Sheet mode in MCS 4 except for the GenerateFASTQ workflow. If using an incompatible sample sheet, import the sample sheet and modify the content in Local Run Manager prior to saving the run. LRM v3-compatible sample sheet templates are available for download from the Illumina support site.

- ▶ **[SupportedModules]**—In this section, specify analysis modules by name, which associates the new kit with an analysis module on the Create Run page. The module name must be entered exactly as shown with no spaces.

Module Name	Description
Assembly	Assembly analysis module
AmpliconDSWorkflow	Amplicon DS analysis module
DNAmpliconWorkflow	DNA Amplicon analysis module
DNAEnrichmentWorkflow	DNA Enrichment analysis module
GenerateFastQWorkflow	Generate FASTQ analysis module
LibraryQCWorkflow	Library QC analysis module
MetagenomicsWorkflow	Metagenomics analysis module
PCRAmplicon	PCR Amplicon analysis module
ResequencingWorkflow	Resequencing analysis module
RNAmpliconWorkflow	RNA Amplicon analysis module
RNAFusionWorkflow	RNA Fusion analysis module
SmallRNAWorkflow	Small RNA analysis module

- 4 Save and rename the file. Do not include spaces in the file name.

## Add a Library Prep Kit

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **Library Prep Kits**.
- 2 Select **Add Library Prep Kit**.
- 3 Navigate to the location of the library kit file and add the file.  
The kit name, description, and read type appears on the Library Prep Kits list.


## Delete a Library Prep Kit

- 1 From the dashboard navigation bar, select the Tools drop-down list, and then select **Library Prep Kits**.
- 2 Select **Delete** next to the kit you want to delete.  
Only custom kits can be deleted from Local Run Manager.
- 3 Select **Delete** to confirm the removal of the kit from Local Run Manager.

## Workflow Overview

This section provides an overview of the workflow for creating a run and viewing analysis results in the Local Run Manager software. Specific steps vary for each analysis module. See the user guide for the analysis module you are using for step-by-step instructions.

### Local Run Manager Workflow

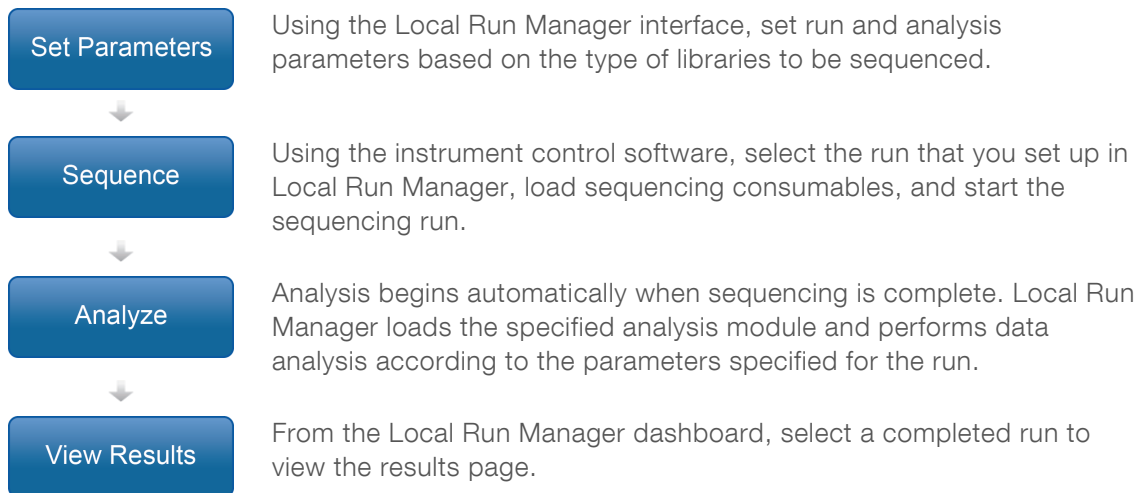

### Set Parameters Overview

To set up a run, select **Create Run** from the Local Run Manager dashboard, and then select an analysis module from the drop-down list. The Create Run page provides options for the selected analysis module. The Create Run page consists of text fields, drop-down lists, toggle settings, and a samples table that is specific to each analysis module. Create Run pages include the following sections:

- ▶ Run Name
- ▶ Run Settings
- ▶ Module-Specific Settings
- ▶ Samples

Fields marked with an asterisk (\*) are required.

### Run Name and Description

The run name is the name that identifies the run from sequencing through analysis. A run name can have up to 40 alphanumeric characters. Spaces and the following special characters, ~!@#\$\$%-\_{}., are allowed. The run name must be unique, and cannot be a name used for a previous run set up. A run description is optional and can have up to 150 alphanumeric characters.

### Run Settings

Run settings identify the following parameters for the run.

- ▶ The index and library prep kits used to prepare samples for sequencing
- ▶ The read type: single read or paired end



- ▶ The number of index reads: 0, 1, or 2
- ▶ The number of cycles for each read
- ▶ Custom primer information (not applicable for off-instrument and iSeq implementations of Local Run Manager)

For some analysis modules, only a default selection is allowed.

For index and library kit compatibility, see the [Local Run Manager support page](#) on the Illumina website.

## Module-Specific Settings

The module-specific settings that appear in this section depend on the analysis module selected for the run. Some modules include options for selecting an aligner and variant caller other than the default methods and configurable analysis settings. Some analysis modules do not include any module-specific settings.

For more information, see the workflow guide for the analysis module you are using.

## Samples Section

For each run, identify samples to be sequenced in the samples table. The samples table includes the sample ID (The Sample ID field has a 40 character limit), an optional sample name, an optional description, plate well (for some index kits), and the indexes used with each sample.

The index columns provide a drop-down list of compatible index sequences based on the library kit specified for the run. For runs using custom index sequences, you can enter the sequence as free text.



### NOTE

During analysis, the iSeq™ 100, MiniSeq, and NextSeq Systems automatically reverse complement the i5 indexes in custom library prep kits. Make sure that the i5 indexes are in the forward orientation.

For analysis modules that require a manifest, the samples table includes a drop-down list of available manifests associated with the module. You can upload and associate a manifest for an analysis module using the Modules & Manifests menu. See [Add Manifest Files to an Analysis Module on page 19](#). To import a manifest for a specific run, use the Import Manifests command.

For analysis modules that perform alignment or other methods that require a reference genome, the samples table includes a drop-down list of available genome folders. See [Set up a Custom Reference Genome on page 26](#).

You can import samples from a compatible text file and export samples from the samples table for use with another run.

Since sample sheet information can vary on a module-to-module basis, see the workflow user guide for the analysis module you are using for module-specific sample sheet settings.

## Import and Export Sample Sheets

If you already have a sample sheet that you want to use to configure the run, select **Import Sample Sheet** to select the CSV file that contains all the run settings and sample information. Make sure that the sample sheet file that you are importing has the same column headings as the sample table for the current analysis module. For example, if the analysis module requires a manifest file, the sample sheet file must have a manifest column.

After filling out the run information and entering at least one sample in the samples table, you can export the run information to an external CSV file by selecting **Export Sample Sheet**. This is helpful to ensure compatibility on future runs or to use as a template when manually editing sample sheets.

**NOTE**

Make sure that you refer to the *Illumina Adapter Sequences document (document # 1000000002694)* and the appropriate analysis module user guide when setting up your sample sheets. Entering the incorrect sequences for your library prep kit will cause your analysis to fail.

## Import Manifest Files for the Run

Unlike importing manifest files on the Modules & Manifests page, importing manifest files on the Create Run imports the manifest for the current run only. Make sure that the manifests you want to import for the run are available in an accessible network location or on a USB drive.

## Available Genomes

In addition to the preinstalled genomes, genome sequence files and reference annotation for other commonly used model organisms are available from the [Illumina iGenomes page](#).

The sequence and annotation files for each iGenome are provided in a compressed file format, \*.tar.gz. Make sure that hg19 is the reference genome when using annotation files. Refer to the iGenomes Overview for installation instructions.

## Set up a Custom Reference Genome

In order to use a custom reference genome in Local Run Manager, the genome files must be set up with a specific folder structure.

- 1 In C:\Illumina\Genomes, create a folder named for the species of your reference genome.
- 2 In the species folder, create a folder named for the agency that published the reference genome. For example, UCSC, NCBI, or Illumina.
- 3 In the agency folder, create a folder named for the version of the reference genome. For example, hg19, mm9, or 2005-10-07.
- 4 In the version folder, create a folder named WholeGenomeFasta.
- 5 Rename the reference genome FASTA file to genome.fa.
- 6 Add genome.fa and its associated GenomeSize.xml file to the WholeGenomeFasta folder.  
The path to the reference genome is:  
C:\Illumina\Genomes\

## View Analysis Results Overview

To view analysis results, select the run name from the Local Run Manager dashboard. A summary of analysis results is presented on the following tabs:

- ▶ Run Overview
- ▶ Sequencing Information
- ▶ Samples & Results

## Run Overview Tab

The Run Overview tab lists information about the run, a summary of sequencing metrics, and the location of the run folder.

Section Title	Description
Run Name / Run ID	The run name assigned when the run was created.
Created By	The name of the user who created the run.
Description	The run description, if provided.
Output Run Folder	The path to sequencing run output folder. Select <b>Copy to Clipboard</b> for quick access to the output folder.
Total Clusters	The number of clusters in a run.
% Clusters PF	The percentage of clusters passing filter.
% ≥ Q30 (Read 1)	The percentage of reads in Read 1 with a quality score of 30 (Q30) or greater.
% ≥ Q30 (Read 2)	The percentage of reads in Read 2 with a quality score of 30 (Q30) or greater.
Last Scored Cycle	The last cycle in the run with a quality score. Unless the run ended early, this value represents the last cycle in the run.

## Sequencing Information Tab

The Sequencing Information tab provides a summary of sequencing run parameters. The following information is provided on the Sequencing Information tab.

Section Title	Description
Instrument Name	The name of the instrument on which the run was performed.
Module Version	The version of the analysis module assigned to the run.
Sequenced By/Imported By	The name of the user who started or imported the run.
Sequencing Start/Imported On	The date and time that the sequencing run began or was imported.
Read Lengths	The number of reads and cycles for each read performed during the run.
Flow Cell Information	The barcode, part number, lot number, and expiration date of the flow cell used for the run.
Buffer Information	The barcode, part number, lot number, and expiration date of the buffer cartridge used for the run.
PR2 Information	The barcode, part number, lot number, and expiration date of the PR2 bottle used for the run.
Reagent Cartridge Information	The barcode, part number, lot number, and expiration date of the reagent cartridge used for the run.

## Samples & Results Tab

The Samples & Results tab provides a summary of analysis results for each sample listed in the samples table, the location of the analysis folder, and an option to requeue analysis.

The following fields are common for all analysis modules. For descriptions of specific analysis results, see the analysis module workflow guide.

Section Title	Description
Select Analysis	A drop-down list of each analysis performed on the data generated from the sequencing run. For each time that analysis is run, a sequential number is assigned. Results are listed separately.
Analysis Folder	Select <b>Copy to Clipboard</b> for quick access to the analysis folder.

## Possible Run States

The Status column on the dashboard shows the status of each run listed. The following table describes the possible states of a run and the appearance of the status bar.

Status	Status Bar Color	Status Description
Ready for Sequencing	None	Waiting for sequencing to begin.
Sequencing Running	Blue	Sequencing in progress.
Sequencing Stopped	Red	Sequencing stopped. No activity.
Sequencing Unsuccessful	Red	Sequencing issue. No activity.
Sequencing Completed	Blue	Sequencing completed. Status bar at 50%.
Primary Analysis Unsuccessful	Red	RTA analysis failed. Status bar at 25%.
Primary Analysis Completed	Blue	Sequencing metrics complete. Status bar at 50%.
Analysis Queued	Blue	Waiting for analysis to begin.
Analysis Running	Blue	Analysis in progress.
Stopping Analysis	Red	Request to stop analysis received.
Analysis Stopped	Red	Analysis stopped. No activity.
Analysis Unsuccessful	Red	Analysis issue. No activity.
Analysis Completed	Green	Activity completed. Status bar at 100%.
Analysis Copying	Blue	Analysis is complete. Data is being copied to the output folder.
Stopping Analysis Copy	Red	Analysis is complete. Request to stop copying of analysis data received.
Copy Stopped	Yellow	Analysis is complete. Copying of analysis data was manually stopped.

## Analysis Module Updates

When a new analysis module or new version of an analysis module is available, you can download the installable file from the Illumina support site and install it onto the Local Run Manager system.

### Install or Update Analysis Modules

- 1 Download the installable file for the new or updated analysis module from the Local Run Manager support page on the Illumina website.
- 2 Copy the installable file (\*.msi or \*.exe) to the instrument or to a USB drive.
- 3 From the instrument computer, double-click the installable file.
- 4 If User Management is turned on, enter your Local Run Manager user name and password when prompted.  
You must have admin-level permissions.
- 5 Select **Next** to begin the installation.
- 6 When the installation is complete, select **Finish**.
- 7 Refresh your browser.
- 8 To confirm that the analysis module is available for use, select the Tools drop-down list, and then select **Modules & Manifests** from the Local Run Manager dashboard.

### Uninstall Local Run Manager

From the computer Control Panel, uninstall components in the following order.

- 1 Uninstall the analysis module. If prompted, enter the Admin user name and password.
- 2 Uninstall Local Run Manager as follows.
  - a Double-click **Local Run Manager**.
  - b Select **Options**.
  - c Select **Uninstall All Components**, and then select **OK**.
  - d Select **Uninstall**.
- 3 Manually delete the reference genome files from their folder locations.

## Revision History

<b>Document</b>	<b>Date</b>	<b>Description of Change</b>
Document # 1000000111492 v01	February 2022	Updated service accounts user name and password requirements to reflect the switch of all instruments to Windows 10, Added uninstallation software guidance, and clarified sample sheet requirements.
Document # 1000000111492 v00	January 2021	Initial release.

## Technical Assistance

For technical assistance, contact Illumina Technical Support.

**Website:** [www.illumina.com](http://www.illumina.com)  
**Email:** [techsupport@illumina.com](mailto:techsupport@illumina.com)

### Illumina Customer Support Telephone Numbers

Region	Toll Free	Regional
North America	+1.800.809.4566	
Australia	+1.800.775.688	
Austria	+43 800006249	+43 19286540
Belgium	+32 80077160	+32 34002973
China	400.066.5835	
Denmark	+45 80820183	+45 89871156
Finland	+358 800918363	+358 974790110
France	+33 805102193	+33 170770446
Germany	+49 8001014940	+49 8938035677
Hong Kong, China	800960230	
Ireland	+353 1800936608	+353 016950506
Italy	+39 800985513	+39 236003759
Japan	0800.111.5011	
Netherlands	+31 8000222493	+31 207132960
New Zealand	0800.451.650	
Norway	+47 800 16836	+47 21939693
Singapore	+1.800.579.2745	
South Korea	+82 80 234 5300	
Spain	+34 911899417	+34 800300143
Sweden	+46 850619671	+46 200883979
Switzerland	+41 565800000	+41 800200442
Taiwan, China	00806651752	
United Kingdom	+44 8000126019	+44 2073057197
Other countries	+44.1799.534000	

Safety data sheets (SDSs)—Available on the Illumina website at [support.illumina.com/sds.html](http://support.illumina.com/sds.html).

Product documentation—Available for download from [support.illumina.com](http://support.illumina.com).



Illumina  
5200 Illumina Way  
San Diego, California 92122 U.S.A.  
+1.800.809.ILMN (4566)  
+1.858.202.4566 (outside North America)  
techsupport@illumina.com  
www.illumina.com

**For Research Use Only. Not for use in diagnostic procedures.**

© 2022 Illumina, Inc. All rights reserved.

**illumina**<sup>®</sup>