

CYP2D6 CNV Caller Software Release Notes

For Infinium Global Screening Array v3

Introduction

These Release Notes detail the known issues present in this initial release of the CYP2D6 CNV Caller software.

I. CYP2D6 CNV Caller

NEW FEATURES:

- N/A

DEFECT REPAIRS:

- N/A

KNOWN ISSUES:

- On Mac OSX, running multiple instances of the CYP2D6 CNV Caller in parallel will error out.
- The software accepts a gencall score cutoff of 0 and completes the analysis when it should throw an error indicating that this is not a valid input.
- The software completes analysis instead of throwing an error when mismatched BPM and CSV manifests are provided as input.
- The software completes analysis instead of throwing an error if a non-csv files is provided for the CSV manifest as input.